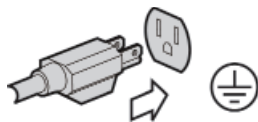
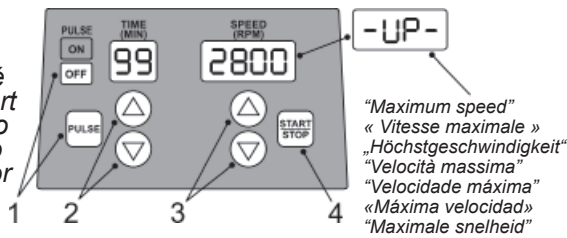
1.0

2.1

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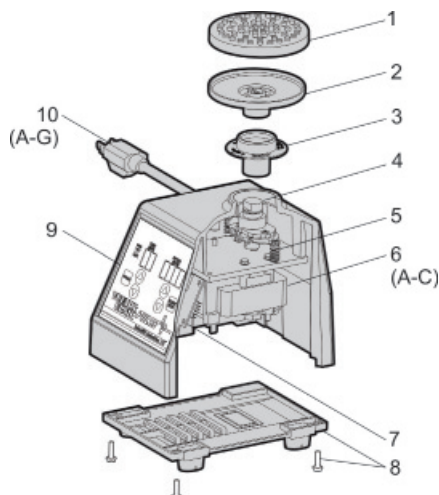
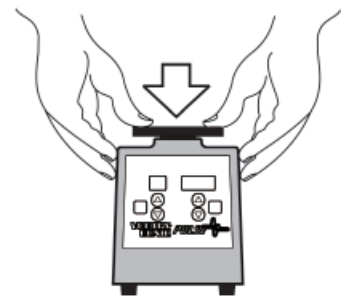
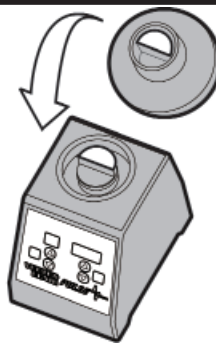
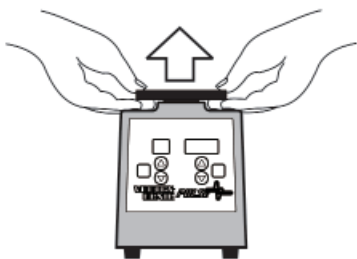
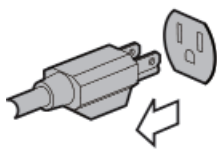
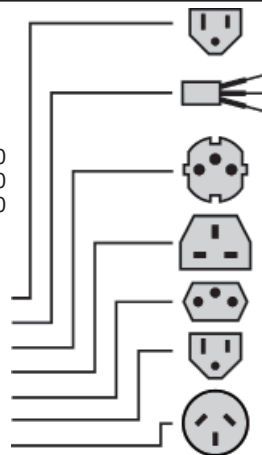
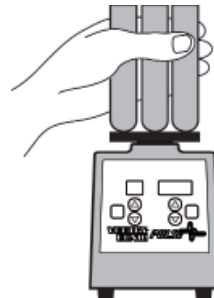
English

Operating Instructions

Retain original packaging for 90 days in case of return. When using the PULSE mode with "On" times of less than 5 seconds, SET speed will be displayed when the unit is mixing, instead of actual speed. Speed control can be adjusted up or down during the mixing cycle, however, TIME can only be adjusted before cycle begins. With prolonged use, your VORTEX-GENIE® PULSE may become hot. This is normal. All VORTEX-GENIE 2 accessories can be used with the VORTEX-GENIE PULSE. *Review instruction 3.0 before attempting to change attachments.*

Caution

- ⚠ If this equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.
- ⚠ Unplug from power before cleaning. Do not immerse.
- ⚠ Do not position the equipment so that it is difficult to disconnect the power cord.
- ⚠ This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

Care and Handling

Your VORTEX-GENIE PULSE should be given the care normally required for any electrical appliance. Unplug unit and wipe housing with damp cloth and detergent. Blot any spills immediately.

Specifications

The VORTEX-GENIE PULSE Mixer is classified as "Installation Category 2", "Overvoltage Category 2" and "Pollution Degree 2".

Environmental: 0°C - 38°C (32°F-100°F), 95% Humidity max. Indoor use only in elevation up to 2000m of altitude.

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